



BELLA BORA

STILL AIR BOX

NEW!

✓ MYCOLOGY

✓ CULTURE
EQUIPMENT

BELLA BORA creates convenient solutions for at-home mycology, combining form and function with superior quality and accessible cost.

DIMENSIONS

THE BELLA BORA STILL AIR BOX IS
2FT LONG X 3FT WIDE X 2 FT HIGH



WHAT IS THE BELLA BORA STILL AIR BOX?

The **BELLA BORA STILL AIR BOX** creates a clean, contaminant-free space for mycology culture work with ample working room to increase the success rate of at home mycologists doing agar transfers, grain inoculation, and grain-to-substrate transfers. The patented design is made with high-quality, durable materials that are scratch-resistant, waterproof, and bleach-safe for easy cleaning. High-visibility plastic and lightweight materials make it ideal for home use.



WHY USE THE BELLA BORA S.A.B.?

The Bella Bora Still Air Box is an essential and effective piece of equipment for mushroom teks involving mushroom spores, liquid culture, agar plates, and sterile inoculation. For the home cultivator or hobbyist interested in taking the next step in mycology, the Still Air Box is significantly more affordable and takes up much less space than a laminar flow hood.

It is superior to the DIY plastic bin SABs, offering a larger workspace at 2ft L x 3ft W x 2 ft H, a convenient arm port design for better mobility and comfort, and side-loading zippers for easy prep. When you're finished, the Bella Bora Still Air Box can be neatly stored in the included case.



QUICK SETUP & PACKS
DOWN FOR TRANSPORT




HOW TO USE BELLA BORA S.A.B.

1. Assemble the still air box according to the included instructions.
2. Get your gloves on, roll up your sleeves, and spray down hands and arms with 70% isopropyl alcohol spray.
3. Prepare your Still Air Box by spraying and wiping down with isopropyl alcohol 70% spray and allow the box to sit while closed for at least 5 minutes.

NOTE: If using isopropyl alcohol, do not utilize open flame in or around the area of the Bella Bora.


You can also use a 10% Bleach solution spray as an alternative to isopropyl alcohol without damaging the Bella Bora.

4. Spray and wipe any items you are using for your clean work before placing them into the still air box.
5. Close the box, spray your gloves and forearms again, and begin your clean work!

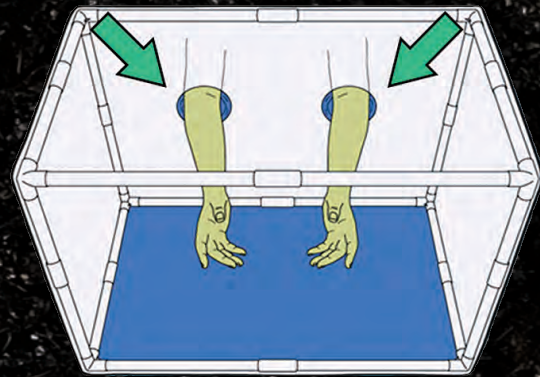

ELASTIC ARM PORTS

Flexible seal allows for a wide range of motion while maintaining a closed environment

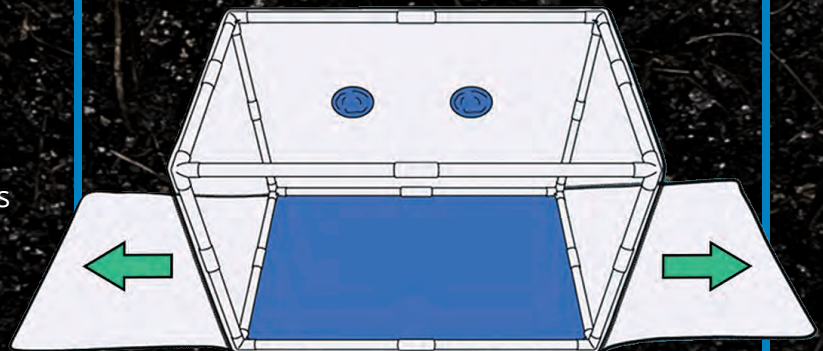

DOUBLE SIDE DOORS

Large zipper opening for convenient loading



COLLAPSIBLE DESIGN

5 minute assembly and disassembly for easy storage and portability